# FINO BALANCE MINI Taschenwaage

REF 51338

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**REF** 51338

**ENG** Instructions for use

 Please read carefully before use  
Keep for referral purposes

**1. Fields of application**  
Compact portable pocket scale for precise weighing for all common weighing units.

**2. Safety precautions**  
No particular hazards worth mentioning.

**3. Processing instructions**

 Information regarding operation:  
Never load the scale beyond its maximum capacity. Overloading can cause permanent damage. Place the objects to be weighed gently on the weighing table. Avoid shaking the scale or letting it fall or any other shocks that could permanently damage the internal sensors. The scale is a precision instrument and must be treated with utmost care.

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Take the scale out of the included pouch and place it on an even, horizontal and clean surface.

Before use leave the scale for a minimum of one hour to adapt to the ambient temperature. Firstly the scale should be calibrated. Refer to section 3.5. Allow the scale sufficient start-up time usually 30 seconds to stabilise the internal components before calibration.

**3.2 Taking measurements**  
Initially the scale is switched on by using the „ON / OFF“- key.

Wait until "0.00" shows in the display. This indicates that the scale is stabilised.

Now place the object to be weighed on the weighing table. The weight that was established by the scale is displayed.

Using the [UNITS]-key you can switch between weighing units.

Using the [LIGHT]-key the background lighting can be switched on or off.

**3.3 Weighing with tare function**  
Firstly the scale is switched on as described in section 3.2. The object to be tared, e.g. a container is placed on the weighing table and the „TARE“-key is pushed. The display switches back to „0.00“. Now the object to be weighed is added and only this weight is displayed as wanted.

**3.4 Adding of weights**  
Switch the scale on and wait until the LCD shows "0.00".

Place a sample quantity (25 pcs. 50 pcs. 75 pcs or 100 pcs.) of the quantity that you want to count on the weighing table.

Press and HOLD the [PCS]-key for 2 seconds. Now the display shows a random number, e.g. "100".

Press the [UNITS]-key, to select the sample quantity on the weighing table.

For confirmation press the [PCS]-key. Now "pcs" is shown in the corner of the display. The scale is now in counting mode.

Remove the sample quantity from the weighing table.

Now you can place any quantity on the scale. The number is shown on the display.

With [UNITS]- or [PCS]-key you can switch from weighing mode to counting mode.

**3.5 Calibration**  
The scale comes calibrated from the manufacturer. Nevertheless the scale should be calibrated regularly and before first use. The used check weight should be calibrated if possible. The weighing unit must be set to grams.

Switch the scale on using the [ON/OFF]-key.

Wait until the LCD shows "0.00" and press now the [UNITS]-key for 3 seconds until the display shows "CAL". Now release the [UNITS]-key.

Press the [UNITS]-key again. The LCD displays "CAL" and subsequently the required check weight.

Place the required check weight. Now the display shows the next check weight (200 g). Add another 100 g check weight or replace the 100g weight with a 200 g check weight.

The display shows "PASS" and switches back to normal weighing mode.

Remove the check weight and press [ON/OFF] to switch the scale off.

The value is stored and the calibration is terminated!

Check the scale by weighing reference weights. It might happen that calibration was not successful. In this case calibration must be repeated.

The variation of 0.01 g might be caused by external factors such as air draughts.

**4. Storage**  
The unit should be stored in a dry place protected from dust and vibrations. Draughts and the direct vicinity to other electronic equipment might negatively influence the precision and reliability of the scale. Avoid extreme heat or cold. If you store the unit for a longer time please remove batteries.

**5. Physical data**  
External dimensions: 11.5 x 6.0 x 2.0 cm; Weighing capacity: 200 g; Readability: 0.01 g; Weighing units g/oz/ozt/dwt; Tare range: Tare fully loaded; Automatic switch-off: after 30 seconds; Operational temperature: 10-30° C; Power supply: 2 x batteries AAA.

**6. Delivery forms**  
FINO BALANCE MINI pocket scale  
each 51338

**7. Guarantee**  
Our technical recommendations of application are based on our own experiences and tests and should only be regarded as guidelines. It rests with the skills and experience of the user to verify that the products supplied by us are suitable for the intended procedures. Our products are undergoing a continuous further development. We reserve the right of changes in construction and composition. It is understood that we guarantee the impeccable quality of our products.

**8. Disposal in accordance with EU guideline 2002/96/EU**

 This equipment may not be disposed in domestic waste! Please contact the manufacturer or your dealer!

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