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9.1 Registrierungsnummer

WEEE-Reg.-Nr. DE 75542610

GUARANTEED TOP QUALITY THROUGHOUT

English

User Information

FINO Stereo Microscopes
86010

By purchasing this product you have made an excellent choice.

FINO quality products from DT&SHOP convince through consistently high quality for economical prices.

1. FIELDS OF APPLICATION

FINO Stereo Microscopes are used for the surface examination in different fields of dental technology.

2. ADVANTAGES

- ◆ 45° slant view
- ◆ Adjustable eye distance
- ◆ Fatigue-free working

3. PHYSICAL DATA

Voltage: 230 V 50 Hz
External dimensions: 19x 11,5x 33
Broad field pair of oculars WF 10x,
Rotating revolver 1x and 3x for
magnification 10x / 30x
Power cable: 1,5 m

4. PROCESSING / INSTRUCTIONS

4.1 Getting started

Please check the unit for transportation damages immediately after unpacking. All transportation damages have to be claimed promptly.

After checking that the voltage on the type label of the unit corresponds with the available mains supply the plug of the unit is connected to the power socket. Place the black rubber ocular shells on the ocular.

4.2 Operation

4.2.1 General operation

To adjust the eye distance both oculars can be swung. The height of the microscope head can be adjusted with black turning knobs to obtain a sharp display of the object to be examined. After loosening the silver locking screws the microscope head can be freely turned. The object magnification (1x or 3x) can be selected by turning of the rotating revolver. Illumination is switched on with the left toggle switch. With the right toggle switch you can select whether the object is illuminated from above or from below. The included object tweezers can be mounted to one of the holding clamps if desired.

4.2.2 Changing of the object tray

Push the tension clamp from the object tray. In the backward area of the microscope is a catch. The object tray is easy to remove. On the right side of the microscope is a small screw, which is used to lock the object tray for transport purposes.

The spring-loaded hinged arm is adjusted to the desired position to ensure relaxed working. Now the lens magnification 1x or 2x can be selected. The eye distance is individually adjustable. If required the top-down illumination can be switched on and can be adjusted to the object.

4.3 Maintenance

To replace the upper bulb remove the silver screw on the left side of the light. The light cover comes loose. The bulb can simply be removed.

To replace the lower bulb remove the glass cover and pull the bulb from the socket.

Exchange the defect bulb with a new 12V / 10W bulb.

On the rear side of the microscope are two fuses. Open the fuse holder with a screwdriver. Replace the defect fuse with a new 250V / 0,5A fuse.

FINO Stereo Microscopes are maintenance-free. Repairs of the unit should only be done by an authorised workshop. The user has to stand good for damages caused by unauthorized/ arbitrary repairs. The unrestricted conditions of delivery and guarantee of FINO apply for this unit.

5. STORAGE

The FINO Stereo Microscopes should be installed in a potential dust-free area and should be covered after use.

6. SAFETY PRECAUTIONS

Do never operate the microscope with wet or moist hands.

7. AVAILABILITY

FINO Stereo-Microscope 86010

With object tweezers, glare-free dark field illumination, dust protection cover, rubber ocular sleeves, spare bulb, spare fuse, object support light / dark.

8. GUARANTEE

Our technical recommendations of application are based on our own experiences and tests and should only be regarded as guidelines. It rests with the skills and experience of the user to verify that the products supplied by us are suitable for the intended procedures. Our products are undergoing a continuous further development. We reserve the right of changes in construction and composition. It is understood that we guarantee the impeccable quality of our products.

9. DISPOSAL IN ACCORDANCE WITH EC GUIDELINE 2002/96/EG

This equipment may not be disposed in domestic waste! Please contact the manufacturer or your dealer!

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